



NEW PIPETMAN® L x1200 MULTICHANNEL MODELS

PIPETMAN® L x1200 models are the latest mechanical multichannel pipettes, combining ergonomics, performance, and secure pipetting.

They cover an extended range of volumes, from 100 to 1200 μL , in either 8- or 12-channel models, helping you increase your daily productivity without sacrificing comfort.

PIPETMAN L x1200 multichannel models are fully autoclavable without disassembly for maximum convenience and protection from cross-contamination.



This leaflet details the specific information related to the new PIPETMAN L x1200 μL multichannel models, in 8- and 12-channels. Refer to the PIPETMAN L user's guide LT801575 for more information.

NOTE

After a maintenance or decontamination procedure, perform a leak test. Refer to page 13 in the PIPETMAN L user's guide.

MODEL	P/N	PIPETMAN DIAMOND TIPS	VOL. (μL)	MAXIMUM PERMISSIBLE ERRORS			
				Gilson		ISO 8655	
				Systematic error (μL)	Random error (μL)	Systematic error (μL)	Random error (μL)
P8x1200L*	FA10039	D1200 DF1200ST	100	± 6.00	≤ 0.9	± 32	≤ 12
			120	± 6.00	≤ 0.9	± 32	≤ 12
			600	± 8.00	≤ 1.2	± 32	≤ 12
P12x1200L**	FA10040		1200	± 10.00	≤ 1.5	± 32	≤ 12

*P8x1200L | 0.235 kg | 265 mm (length)

**P12x1200L | 0.285 kg | 265 mm (length)

SPARE PARTS

LABEL	DESCRIPTION	PART NUMBER 8X1200	PART NUMBER 12X1200
1	Tip ejector	FA07099	FA07100
2	Ejector lock	F507008	F507008
3	Ejector spacer	F507001	F507003



WARRANTY

Gilson warrants this pipette against defects in material under normal use and service for a period of **three years** from the date of purchase. This warranty shall not apply to pipettes which are subject to abnormal use and/or improper or inadequate maintenance (contrary to the recommendations given in the user's guide), including, but not limited to pipettes which have been subjected to physical damage, improper handling, or spillage or exposure to any corrosive environment. This warranty shall also be void in the event pipettes are altered or modified by any party other than Gilson or its designates. Gilson's sole liability under this warranty shall be limited to, at Gilson's sole option, repair, or replacement of any defective components of pipettes or refund of the purchase price paid for such pipettes. **Routine cleaning, control, and recalibration are not covered under the warranty. The replacement of wearing parts such as seals, O-rings, broken pistons assembly, and broken tip holders are not covered under the warranty.**

NOTICE

Yearly routine maintenance is highly recommended to keep your pipette in good condition, ensuring a continued high level of performance.

THE FOREGOING WARRANTY IS EXCLUSIVE AND GILSON HEREBY DISCLAIMS ALL OTHER WARRANTIES, WHETHER EXPRESS OR IMPLIED, INCLUDING ANY WARRANTIES OF MERCHANTABILITY AND ANY WARRANTIES OF FITNESS FOR A PARTICULAR PURPOSE, UNDER NO CIRCUMSTANCES SHALL GILSON BE LIABLE FOR ANY CONSEQUENTIAL, PUNITIVE, INDIRECT OR INCIDENTAL DAMAGES ARISING OUT OF ANY BREACH OF ANY EXPRESS OR IMPLIED WARRANTY.

www.gilson.com/contactus

LT801633/A | ©2019 Gilson, Inc.
All rights reserved.

 **GILSON**[®]